



Step 1.

Place item to be tested on a flat and clean surface.

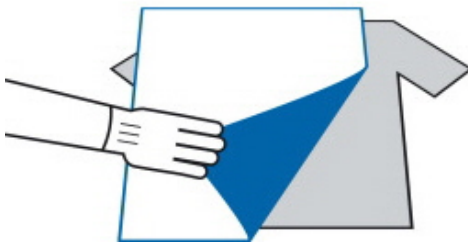


Step 2.

Dampen Phadebas® paper with water. Cover testing area with Phadebas® paper, reagent side down.

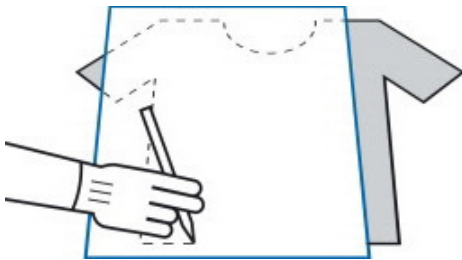


Step 3.

Trace an outline of the testing area on the Phadebas® paper.

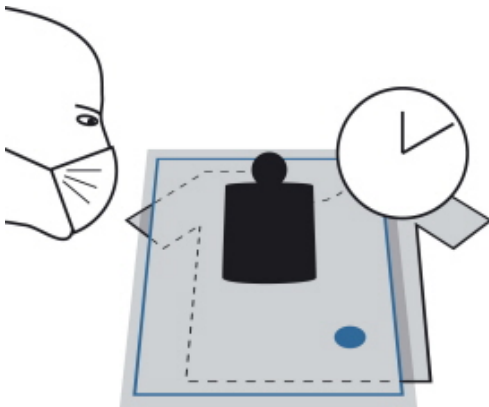

Step 4.

Cover the Phadebas® paper with a clean glass board. Weigh down the arrangement.

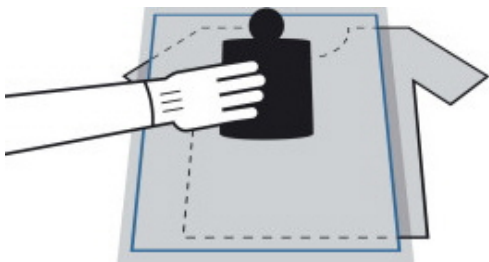

Step 5.

Start the timer and observe the test frequently during 40 minutes. Record the time of positive reaction (distinct area of diffuse blue color).

For a detailed operating procedure, visit www.phadebas.com/archive/