




DETECTS & IDENTIFIES SALIVA STAINS


OPERATING PROCEDURE

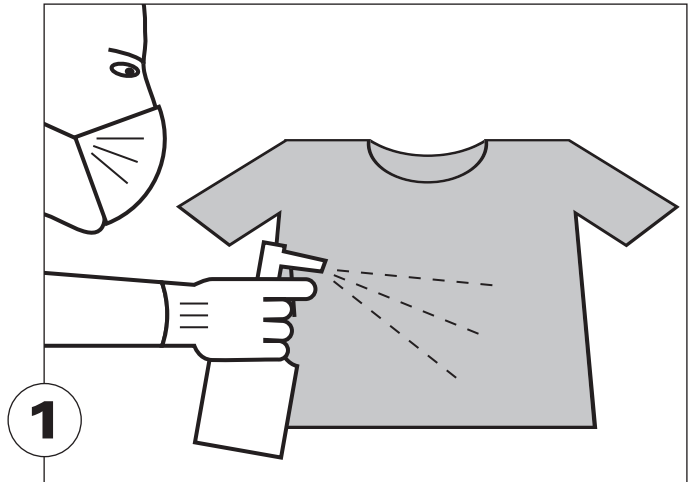


Place item to be tested on a flat and clean surface. Dampen the item with water.




- Store in closed casing at room temperature.
- Do not remove desiccants.
- Avoid contamination – wear exam gloves and mouth mask.

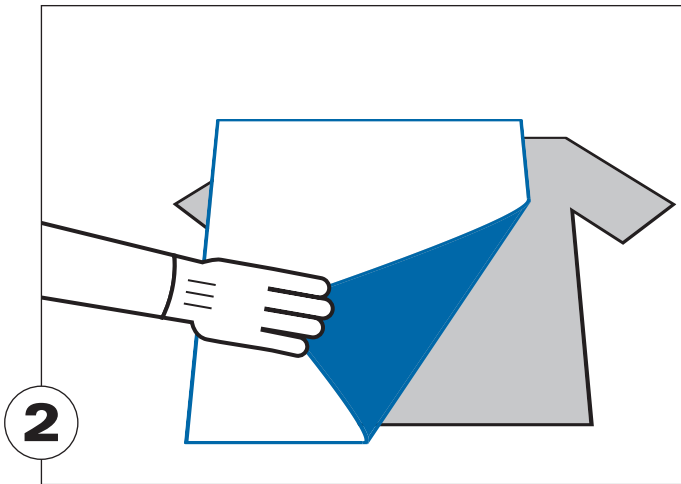


Cover testing area with Phadebas® paper, reagent side down. Dampen Phadebas® paper with water.

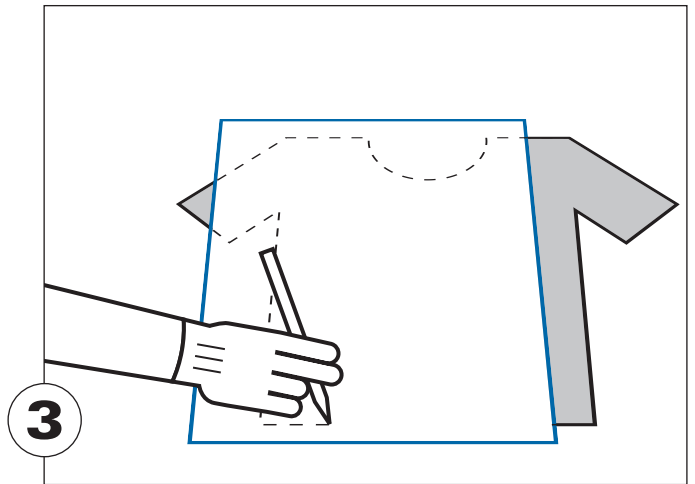

Trace an outline of the testing area on the Phadebas® paper.

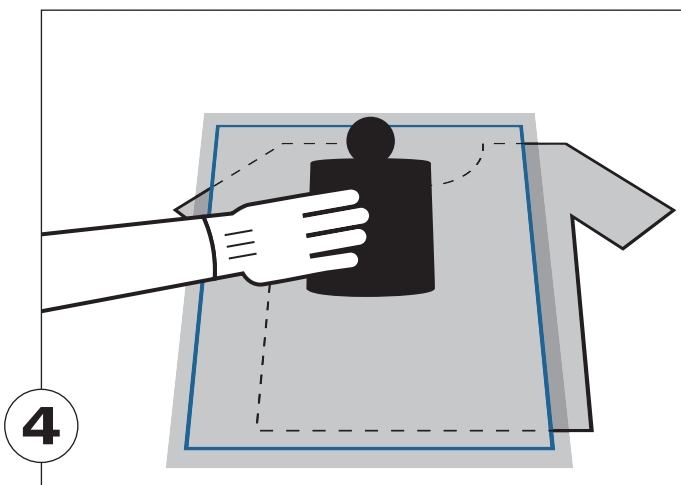

Cover the Phadebas® paper with a clean glass board. Weigh down the arrangement.

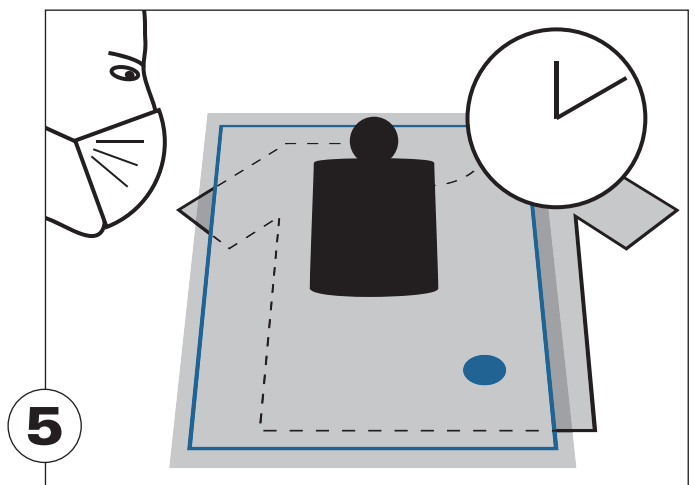

Start the timer and observe the test frequently during 40 minutes. Record the time of positive reaction (distinct area of diffuse blue color).