



**Step 1.**

Place item to be tested on a flat and clean surface.

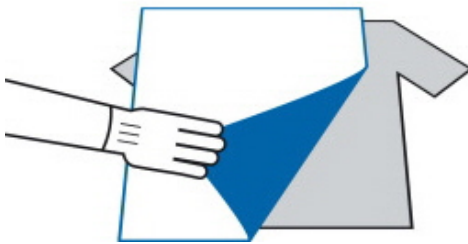


**Step 2.**

Dampen Phadebas® paper with water. Cover testing area with Phadebas® paper, reagent side down.

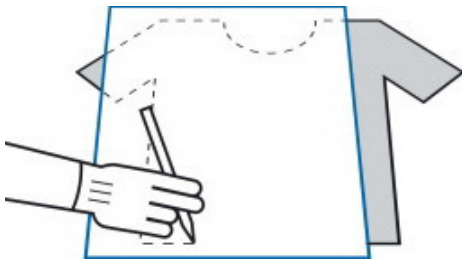


**Step 3.**

Trace an outline of the testing area on the Phadebas® paper.

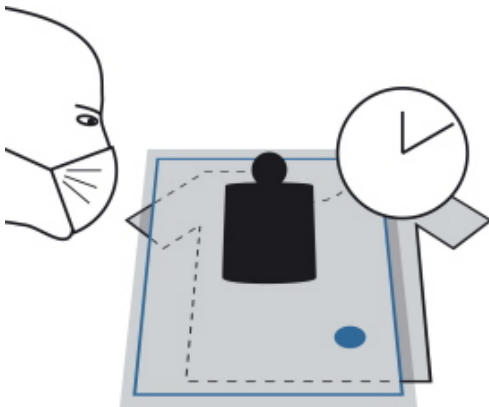

**Step 4.**

Cover the Phadebas® paper with a clean glass board. Weigh down the arrangement.

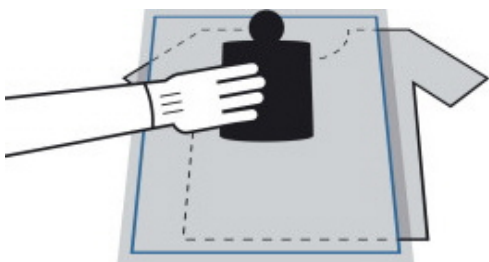

**Step 5.**

Start the timer and observe the test frequently during 40 minutes. Record the time of positive reaction (distinct area of diffuse blue color).

For a detailed operating procedure, visit [www.phadebas.com/archive/](http://www.phadebas.com/archive/)